

PURE-EVs

Size exclusion chromatography columns for extracellular vesicle isolation

Product Code: HBM-PEV-##. Quantity: 5 or 10 SEC columns

PURE-EVs Columns

Size Exclusion Chromatography (SEC) is considered one of the best methods for isolating and purifying Extracellular Vesicles (EVs) from different matrices. This technique is very efficient for separating EVs from the circulating proteins and does not affect the original shape and functionality of the vesicles.

PURE-EVs is a SEC column designed for isolating EVs in a fast and easy way from the volume amounts suggested:

Fluid	Volume amount
Plasma	0.2 ml up to 2 ml
Serum	0.2 ml up to 2 ml
Urine	0.2 ml up to 2 ml *
Cell media	0.2 ml up to 2 ml *

* For diluted fluids concentration 10-fold is recommended.

Procedure for EV isolation

1. Sample preparation.


Prepare the sample by centrifugation steps as suggested in the table below:

Fluid	Suggested	Optional
Plasma	10 min at 300 g (save super) 20 min at 1200 g (save super)	30 min at 10000 g (to eliminate vesicles > 200 nm)
Serum	10 min at 300 g (save super) 20 min at 1200 g (save super)	30 min at 10000 g (to eliminate vesicles > 200 nm)
Urine*	10 min at 300 g (save super).	
Cell media*	10 min at 300 g (save super) 20 min at 1200 g (save super).	30 min at 10000 g (to eliminate vesicles > 200 nm).

* Biofluids which present a diluted population of EVs, can be concentrated 10 folds. Viscous fluids (saliva) must be diluted in PBS 1x before to proceed to the EV isolation in SEC columns. Use of TFF-Easy for concentrating the diluted matrices is recommended. Alternatively MWCO concentrators (100K) can be used.


2- Column preparation.

- PURE-EVs columns are provided with a layer of preservative buffer.
- Open the upper and the lower cap of the PURE-EVs column and let to flow almost all the buffer throught the column, avoiding to dry the surface of the gel.
- Wash the column with 3 volumes of PBS 1x buffer (3 x 10 ml) to eliminate preservative buffer residues.

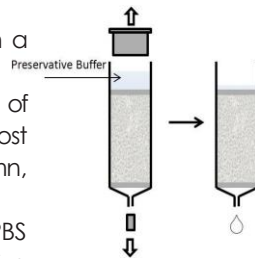


3- Sample loading.

- Rinse the column with 200 µl up to 2 ml of sample containing EVs.
- Collect 500 µl fractions.
- When the sample is inside the gel matrix rinse the column with PBS 1x. PBS 1x is the mobile phase of SEC column, do not let the column to get dried.

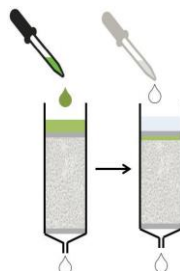

4- EV isolation.

Separation of EV and circulating proteins proceeds as indicated in the figure below (collecting 500 µl fractions).

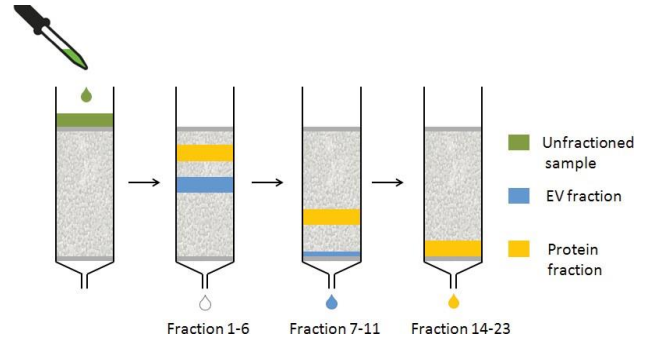


5- Column washing.

- After all fractions are collected wash the column from the residues of sample with approximately 20 ml of PBS. Never get the column dried. After the last washing step add to the column 1 ml of PBS 1x and close the caps. Column can be stored at 4°C and reused up to 5 times.

Results and EV Separation

PURE-EVs column was filled with 2 ml of medium from HCT116 cells, concentrated 10 folds by TFF-Easy. 24 fractions (500 µl each one) have been collected and analyzed by ELISA ExoTEST™ assay (CD81 marker) and by BCA test for determining respectively vesicle and total protein content. EVs are eluted in fractions 8 - 12 (turnaround time approximately 15 min), whereas circulating proteins corresponded to the fractions 14 - 24.

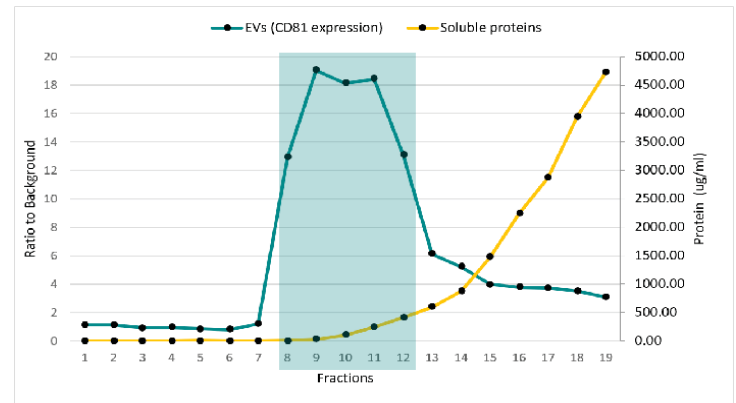

EV Isolation: FAST protocol


The FAST protocol allows to obtain EV preparation in approximately 15 minutes, without collecting all the 500 µl fractions.

- Prepare the samples as indicated in "Sample preparation" paragraph.
- Prepare the column as indicated in "Column preparation" paragraph.
- Fill the column with 200 µl up to 2 ml of sample containing EVs.
- Collect and discard the first 3 ml fraction (void volume), which does not contain vesicles*.

*When the sample is inside the gel matrix rinse the column with PBS 1x. PBS 1x is the mobile phase of SEC column, do not let the column to get dried.




- Collect and save the second 2 ml fraction, containing EVs.
- After EVs are collected wash the column as indicated in "Washing column paragraph". Store the column at 4 °C.

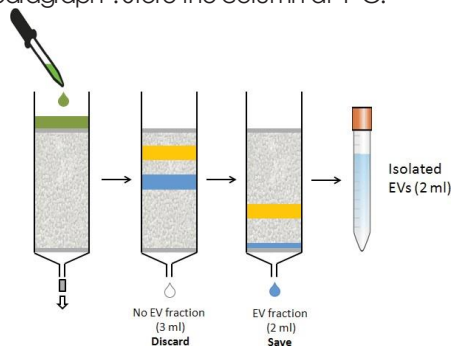

EV Separation with FAST Protocol

Collected fractions were analyzed by NTA (Zetaview, Particle Metrix), ExoTEST and by BCA assay for determining EV and total protein content. EVs are eluted in fraction 2.

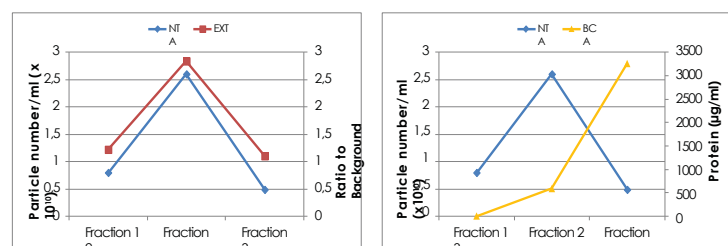

EV elution peak. ExoTEST™ vs NTA analysis.

EV elution vs circulating protein elution. NTA analysis compared to protein BCA test.

References

Apostolov A., Mladenović, D., Tilk, K., Löhmus, A., Vesselin Baev, Yahubyan, G., Sola-Leyva, A., Bergamelli, M., André Görgens, Zhao, C., Samir E L Andaloussi, Aive Kalinina, Acharya, G., Lanner, F., Merli Saare, Peters, M., Piomboni, P., Luddi, A., Salumets, A., & Aleksejeva, E. (2025). Multi-omics analysis of uterine fluid extracellular vesicles reveals a resemblance with endometrial tissue across the menstrual cycle: biological and translational insights. *Human Reproduction Open*, 2025(2). <https://doi.org/10.1093/hropen/hoaf010>

Ikeda, H., Kawase, K., Nishi, T., Watanabe, T., Takenaga, K., Inozume, T., Ishino, T., Aki, S., Lin, J., Kawashima, S., Nagasaki, J., Ueda, Y., Suzuki, S., Makinoshima, H., Itami, M., Nakamura, Y., Tatsumi, Y., Suenaga, Y., Morinaga, T., . . . Togashi, Y. (2025). Immune evasion through mitochondrial transfer in the tumour microenvironment. *Nature*. <https://doi.org/10.1038/s41586-024-08439-0>

Hass, R., Von Der Ohe, J., & Luo, T. (2024). Human mesenchymal stroma/stem-like cell-derived taxol-loaded EVs/exosomes transfer anti-tumor microRNA signatures and express enhanced SDF-1-mediated tumor tropism. *Cell Communication and Signaling*, 22(1). <https://doi.org/10.1186/s12964-024-01886-2>

Kim, H. K., Choi, Y., Kim, K. H., Byun, Y., Kim, T. H., Kim, J. H., An, S. H., Bae, D., Choi, M. K., Lee, M., Kang, G., Chung, J., Kim, S., & Kwon, K. (2024). Scalable production of siRNA-encapsulated extracellular vesicles for the inhibition of KRAS-mutant cancer using acoustic shock waves. *Journal of Extracellular Vesicles*, 13(9). <https://doi.org/10.1002/jev2.12508>

Santoro, J., Nuzzo, S., Franzese, M., Salvatore, M., & Grimaldi, A. M. (2024). Goat milk extracellular vesicles: Separation comparison of natural carriers for theragnostic application. *Heliyon*, 10(6), e27621. <https://doi.org/10.1016/j.heliyon.2024.e27621>

Aparicio, P., Navarrete-Villanueva, D., Gómez-Cabello, A., López-Royo, T., Santamaría, E., Fernández-Irigoyen, J., Ausín, K., Arribeo, M., Sebastian, V., Vicente-Rodríguez, G., Osta, R., & Manzano, R. (2024). Proteomic profiling of human plasma extracellular vesicles identifies PF4 and C1R as novel biomarker in sarcopenia. *Journal of Cachexia Sarcopenia and Muscle*. <https://doi.org/10.1002/jcsm>

Hamdi, M., Sánchez, J. M., Fernandez-Fuertes, B., Câmara, D. R., Bollwein, H., Rizo, D., Bauersachs, S., & Almiñana, C. (2024). Oviductal extracellular vesicles miRNA cargo varies in response to embryos and their quality. *BMC Genomics*, 25(1). <https://doi.org/10.1186/s12864-024-10429-5>

Nebogatova, J., Härk, H. H., Puskar, A., Porosk, L., Guazzi, P., Dowaidar, M., Langel, Ü., & Kurikoff, K. (2023). A Method for Using Cell-Penetrating Peptides for Loading Plasmid DNA into Secreted Extracellular Vesicles. *Biomolecules*, 13(12), 1751. <https://doi.org/10.3390/biom13121751>

Liu, C., Seneviratne, C. J., Palma, C., Rice, G., Salomon, C., Khanabdal, R., Ivanovski, S., & Han, P. (2023). Immunoaffinity-enriched salivary small extracellular vesicles in periodontitis. *Extracellular Vesicles and Circulating Nucleic Acids*, 4(4), 698–712. <https://doi.org/10.20517/evcna.2023.48>

Misawa, T., Hitomi, K., Miyata, K., Tanaka, Y., Fujii, R., Chiba, M., ... & Takahashi, A. (2023). Identification of Novel Senescent Markers in Small Extracellular Vesicles. *International journal of molecular sciences*, 24(3), 2421.

Casado-Díaz, A., Jiménez-Navarro, C., Torrecillas-Baena, B., Camacho-Cardenosa, M., Quesada-Gómez, J. M., & Gálvez-Moreno, M. Á. (2023). Effect of extracellular vesicles derived from hypoxia-preconditioned human mesenchymal stem cells on osteoblastogenesis and adipogenesis in vitro. *Revista De Osteoporosis Y Metabolismo Mineral*. <https://doi.org/10.20960/revosteoporosmetabminer.00012>

Sakr, O. G., Gad, A., Cañón-Beltrán, K., Cajas, Y. N., Prochazka, R., Rizo, D., & Rebollar, P. G. (2023). Characterization and identification of extracellular vesicles-coupled miRNA profiles in seminal plasma of fertile and subfertile rabbit bucks. *Theriogenology*, 209, 76–88. <https://doi.org/10.1016/j.theriogenology.2023.06.020>

Hayasaka, R., Tabata, S., Hasebe, M., Ikeda, S., Hikita, T., Oneyama, C., Yoshitake, J., Onoshima, D., Takahashi, K., Shibata, T., Uchida, K., Baba, Y., Soga, T., Tomita, M., & Hirayama, A. (2023). Metabolomics of small extracellular vesicles derived from isocitrate dehydrogenase 1-mutant HCT116 cells collected by semi-automated size exclusion chromatography. *Frontiers in Molecular Biosciences*, 9. <https://doi.org/10.3389/fmolb.2022.1049402>

Lin, S. C., Li, W. N., Lin, S. C., Hou, H. T., Tsai, Y. C., Lin, T. C., ... & Tsai, S. J. (2023). Targeting YAP1 ameliorates progesterone resistance in endometriosis. *Human Reproduction*, 38(6), 1124–1134.

Zabegina, L., Zyatchin, I., Kniazeva, M., Shalaev, A., Berkut, M., Sharoyko, V., ... & Malek, A. (2023). Diagnosis of Prostate Cancer through the Multi-Ligand Binding of Prostate-Derived Extracellular Vesicles and miRNA Analysis. *Life*, 13(4), 885.

Jiménez-Navarro, C., Torrecillas-Baena, B., Camacho-Cardenosa, M., Quesada-Gómez, J. M., Gálvez-Moreno, M. Á., & Casado-Díaz, A. Effect of extracellular vesicles derived from hypoxia-preconditioned human mesenchymal stem cells on osteoblastogenesis and adipogenesis in vitro.

Sakr, O. G., Cañón-Beltrán, K., Cajas, Y. N., Rizo, D., & Rebollar, P. G. (2022). Characterization of extracellular vesicles in seminal plasma of fertile and subfertile rabbit bucks. *Animal Reproduction (AR)*, 19(2), 0-0.

Yim, K. H. W., Borgoni, S., & Chahwan, R. (2022). Serum extracellular vesicles profiling is associated with COVID-19 progression and immune responses. *Journal of Extracellular Biology*, 1(4), e37.

Leal, C. L. V., Cañón-Beltrán, K., Cajas, Y. N., Hamdi, M., Yaryes, A., de la Blanca, M. G. M., ... & Rizo, D. (2022). Extracellular vesicles from oviductal and uterine fluids supplementation in sequential in vitro culture improves bovine embryo quality.

Sun, C. Y., Chen, G. D., He, B. C., Fu, W. E., Lee, C. H., Leu, Y. W., & Hsiao, S. H. (2022). Dysregulated HIC1 and RassF1A expression in vitro alters the cell cytoskeleton and exosomal Piwi-interacting RNA. *Biochemical and Biophysical Research Communications*.

Rana, N., Suliman, S., Al-Sharabi, N., & Mustafa, K. (2022). Extracellular Vesicles Derived from Primed Mesenchymal Stromal Cells Loaded on Biphasic Calcium Phosphate Biomaterial Exhibit Enhanced Macrophage Polarization. *Cells*, 11(3), 470.

Nikiforova, N., Chumachenko, M., Nazarova, I., Zabegina, L., Slyusarenko, M., Sidina, E., & Malek, A. (2021). CM-Dil Staining and SEC of Plasma as an Approach to Increase Sensitivity of Extracellular Nanovesicles Quantification by Bead-Assisted Flow Cytometry. *Membranes*, 11(7), 526.

